

**foodproof® *Enterobacteriaceae* plus *Cronobacter* Detection Kit
 - Hybridization Probes (LC 2.0, 480) -**

**Order No. R 310 15.1
 Order No. R 310 15.1 L**

**Quick Reference Procedure (for LC 2.0)
 Version 2, April 2016**

A. LightCycler® Carousel-Based System Protocol

The following procedure is optimized for use with the LightCycler® 2.0 Carousel-Based System. Program the LightCycler® Carousel-Based System before preparing the reaction mixes. Use the following LightCycler® 2.0 System PCR-program for the **foodproof *Enterobacteriaceae* plus *Cronobacter* Detection Kit** (for details on how to program the experimental protocol, see the LightCycler® 2.0 Instrument Operator's Manual):

Pre-incubation (Prevention of carry-over contamination, activation of Taq DNA polymerase DNA-free, denaturation of template DNA)				
Programs/Cycle Program Data		Value		
Cycles	1			
Analysis Mode	None			
Temperature Targets		Segment 1	Segment 2	
Target/Target Temperature [°C]	37	95		
Hold/Incubation Time [h:min:s]	00:04:00	00:02:00		
Ramp Rate/Temperature Transition Rate [°C/s]	20	20		
Sec Target/Secondary Target Temperature [°C]	0	0		
Step Size [°C]	0.0	0.0		
Step Delay [cycles]	0	0		
Acquisition Mode	None	None		
Amplification (of the target DNA)				
Programs/Cycle Program Data		Value		
Cycles	34			
Analysis Mode	Quantification			
Temperature Targets		Segment 1	Segment 2	Segment 3
Target/Target Temperature [°C]	95	64	72	
Hold/Incubation Time [h:min:s]	00:00:02	00:00:35	00:00:20	
Ramp Rate/Temperature Transition Rate [°C/s]	20	20	20	
Sec Target/Secondary Target Temperature [°C]	0	60	0	
Step Size [°C]	0.0	0.2	0.0	
Step Delay [cycles]	0	8	0	
Acquisition Mode	None	Single	None	
Cooling (the rotor and thermal chamber)				
Programs/Cycle Program Data		Value		
Cycles	1			
Analysis Mode	None			
Temperature Targets		Segment 1		
Target/Target Temperature [°C]	40			
Hold/Incubation Time [h:min:s]	00:00:30			
Ramp Rate/Temperature Transition Rate [°C/s]	20			
Sec Target/Secondary Target Temperature [°C]	0			
Step Size [°C]	0.0			
Step Delay [cycles]	0			
Acquisition Mode	None			

Continued next page

B. Preparation of the PCR Mix

Proceed as described below to prepare a 20 µl standard reaction.

Do not touch the surface of the capillaries. Always wear gloves when handling the capillaries.

1. Depending on the total number of reactions, place the required number of LightCycler® Capillaries in centrifuge adapters or in a LightCycler® Sample Carousel in a LC Carousel Centrifuge Bucket.
2. Thaw the solutions and, for maximal recovery of contents, briefly spin vials in a microcentrifuge before opening. Mix carefully but thoroughly by pipetting up and down.
3. In a 1.5 ml reaction tube, prepare the PCR Mix by adding the following components in the order mentioned below, then mix gently but thoroughly by pipetting up and down:

The volumes indicated below are based on a single 20 µl standard reaction. Prepare the PCR mix by multiplying the amount in the "Volume" column by the number of reactions to be cycled plus one or two additional reactions to cover pipetting losses.

Component	Volume
foodproof <i>Enterobacteriaceae</i> plus <i>Cronobacter</i> Master Mix, (vial 1, yellow cap)	13 µl
foodproof <i>Enterobacteriaceae</i> plus <i>Cronobacter</i> Enzyme Solution, (vial 2, red cap)	1 µl
foodproof <i>Enterobacteriaceae</i> plus <i>Cronobacter</i> Internal Control - LC 2.0, (vial 3, white cap)	1 µl
Total volume	15 µl

4. • Mix carefully but thoroughly by pipetting up and down. Do not vortex.
 - Pipet 15 µl PCR mix into each LightCycler® capillary.
 - For the samples of interest, add 5 µl sample DNA to a capillary, seal with a stopper.
 - For the negative control, add 5 µl H₂O PCR-grade (vial 5, colorless cap), seal with a stopper.
 - For the positive control, add 5 µl foodproof *Enterobacteriaceae* plus *Cronobacter* Control Template (vial 4, purple cap), seal with a stopper.
5. • Place the adapters (containing the capillaries) in a standard benchtop microcentrifuge. (Place the centrifuge adapters in a balanced arrangement within the centrifuge.)
 - Centrifuge at 700 x g for 5 s (3,000 rpm in a standard benchtop microcentrifuge).
 - Alternatively, use the LC Carousel Centrifuge for spinning the capillaries.
6. Transfer the capillaries to the LightCycler® Carousel Based System.
7. Cycle the samples as described above.

Note: For further information please refer to: www.bc-diagnostics.com